



Product Description

The CompuFlow 1000 flow pump system is designed to generate realistic, accurate and repeatable physiological volume flow waveforms, including pre-programmed; carotid, femoral, sine, square and constant flow waveforms. The system includes an embedded industrial grade motherboard and proprietary *SimuFlow III Waveform Editing Software* for creating user-defined physiological flow waveforms. Users have the option of inputting their own waveform data points, reshaping the supplied pre-programmed waveforms (click and drag) or downloading waveforms onto the system.

The CompuFlow 1000 is accurate to within $\pm 1.0\%$ over a flow range of 0.1 to 35 ml/s. Each system is calibrated and accompanied with a *certificate of calibration* specifying the calibration results at relevant flow rates, as well as being provided with a calibration constant to verify the pump systems accuracy.

The CompuFlow 1000 is two-unit design, a Control Unit and a Pump Unit with an internal reservoir.

Combine the CompuFlow 1000 with an appropriate vascular phantom and blood mimicking fluid, and the resulting closed flow loop ensures easy, accurate and reliable evaluation and validation of diagnostic imaging systems and techniques. The system is ideal for flow quantification in Doppler ultrasound, CTA, particle image velocimetry (PIV) and endovascular simulation applications.

Applications

Research and new technology developments requiring accurate and repeatable pulsatile and constant flow waveforms (i.e. flow quantification, Doppler ultrasound, CTA, particle image velocimetry (PIV) and endovascular techniques.

Ability to control waveform shape.

Gated flow studies.

Calibration of clinical imaging systems.

Product Features

REALISTIC PHYSIOLOGICAL FLOW WAVEFORMS
PULSATILE AND CONSTANT FLOW WAVEFORMS
ACCURATE VOLUME FLOW WAVEFORMS
REPRODUCIBLE WAVEFORMS
PROGRAMMABLE USER DEFINED WAVEFORMS
CALIBRATION CERTIFICATE
MULTIMODALITY IDEAL FOR COMPARING IMAGING MODALITIES

Accurate and reproducible physiological flow waveforms including those with reverse flow components such as the femoral waveform. Accurate ($\pm 1.0\%$) physiological waveforms.

SimuFlow III waveform editing software allows for the development of user-defined pulsatile and constant flow waveforms. Users can control the shape of the waveform.

Pre-programmed waveforms include; carotid, femoral, sine, rectangular and constant flow waveforms.

Control options include the use of a remote hand-held LCD display/keypad or alternatively a keyboard, mouse and monitor can be connected to control the embedded computer.

Calibration certificate is provided with each system verifying the systems accuracy. The system can be calibrated and certified traceably to a national standard for regulatory requirements. A re-calibration service is available.





Fluid Flow Specifications

Volume range 0.1 – 35 ml/s
Accuracy of $\pm 1.0\%$
Sine waves up to 12 Hz
Physiological flow waveforms including reverse flow components

Mechanical Specifications

High-torque micro-stepping motor
450 ml pump cylinder
Adjustable over-pressure shut-off (7 to 50 psi)
Self-sealing 1/4 inch inner diameter plastic connectors

Computer Control

On-board dedicated motor controller
Easily upgraded high level control software
Windows XP Embedded Operating System

Interface Options

Hand-held remote control keypad with backlit LCD
User supplied monitor, keyboard and mouse

Physical

Control Unit dimensions: 14.0" (L) x 11.0" (W) x 9.0" (H)
Control Unit Weight: 15.0 lbs. (6.8 kg)
Pump Unit dimensions: 19.6" (L) x 10.0" (W) x 9.25" (H)
Pump Unit Weight: 38 lbs. (17.2 kg)
Reservoir capacity: 1.5 litres
Circulating Fluid Volume: 2 litres

File Format

Waveform data is stored on embedded harddrive.
Files can contain up to 1,000 flow-rate values describing a waveform
User specified time-base ranging from 2 to 50 ms per point
Waveforms can be looped continuously
Variable frequency waveforms can be generated

Electrical

110/220 VAC 50/60 Hz, 4.0 amps @ 120V
EMI shielding
1 RS-232 serial communication port
1 remote controller port
1 ECG TTL trigger output (BNC)
Monitor port
Keyboard port
Mouse port
Two cables to connect the Control & Pump Units

Calibration Options

Standard calibration
Traceable calibration to a National Standard

Accessory Products

U-245 Doppler Tissue Flow Phantom
Carotid Bifurcation Models for Doppler Applications
Custom Doppler Tissue Flow Phantoms
Blood mimicking fluids, BMF-US
Multi+ Variable Waveform Software Option to mimic arrhythmia



S H E L L E Y M E D I C A L I M A G I N G T E C H N O L O G I E S

Sales Office

157 Ashley Crescent, London, Ontario N6E 3P9 Canada
Phone: 1 (519) 690-0874 Fax: 1 (519) 690-0875
Email: bob.gravett@simutec.com Web: www.simutec.com

Service Centre

41 Coldwater Road, Toronto, Ontario M3B 1Y8 Canada
Phone: 1 (416) 447-6471 Fax: 1 (416) 447-9313
Email: service@simutec.com

A D I V I S I O N O F S H E L L E Y A U T O M A T I O N I N C .